


## Datasheet

## OE-200-UV

### Variable Gain Photoreceiver – Fast Optical Power Meter

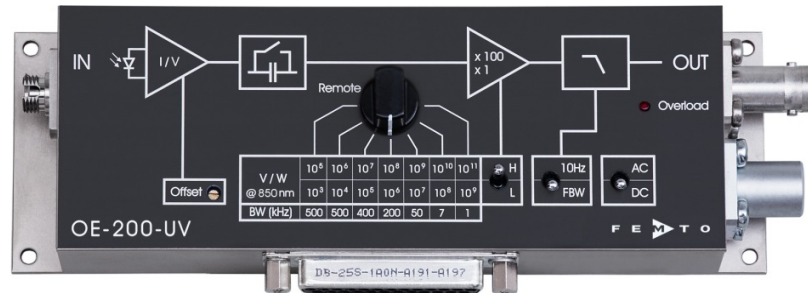

The picture shows model OE-200-UV-FC with fiber optic input.

<p>Features</p>	<ul style="list-style-type: none"> <li>• <b>Si-PIN detector, active area 1.1 x 1.1 mm<sup>2</sup></b></li> <li>• <b>Spectral range 190 - 1000 nm</b></li> <li>• <b>Very low noise, NEP down to 17 fW/√Hz</b></li> <li>• <b>Bandwidth up to 500 kHz</b></li> <li>• <b>Conversion gain adjustable from 1 x 10<sup>3</sup> up to 1 x 10<sup>11</sup> V/W</b></li> <li>• <b>Optical free-space input 1.035"-40 threaded, alternatively 25 mm diameter unthreaded</b></li> <li>• <b>Fiber optic input available as screw-on adapter (1.035"-40) and as permanently mounted FC-input (for calibrated precision measurements)</b></li> <li>• <b>Factory calibrated at 850 nm (fiber optic FC version only)</b></li> <li>• <b>Full manual and remote control capability</b></li> </ul>
<p>Applications</p>	<ul style="list-style-type: none"> <li>• <b>All-purpose very low-noise photoreceiver (O/E converter)</b></li> <li>• <b>Time resolved optical pulse and power measurements</b></li> <li>• <b>Optical front-end for oscilloscopes, spectrum analyzers, A/D converters and lock-in amplifiers</b></li> <li>• <b>Fast fiber optic power meter</b></li> </ul>
<p>Block Diagram</p>	

# Variable Gain Photoreceiver – Fast Optical Power Meter

Available Versions

OE-200-UV-FST




Internal threaded coupler ring with 30 mm outer diameter (included)

1.035"-40 threaded flange for free space applications compatible with many optical standard accessories and for use with various types of fiber connector adapters.

Optional: Fiber adapters PRA-FC and PRA-FSMA



OE-200-UV-FS

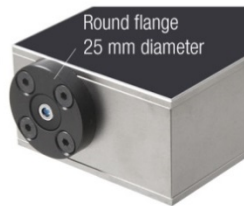

25 mm dia. unthreaded flange for free space applications compatible with many optical standard accessories.

OE-200-UV-FC



fix/permanent FC fiber connector for highest coupling efficiency and best conversion gain accuracy ( $\pm 5\%$ )

Since illumination conditions with the permanently mounted fiber optic connector are well defined, the FC model is delivered with a factory calibrated conversion gain at 850 nm.

The electro optical conversion gain factors of the FST and FS free space models are set to fit nominally at 850 nm.

## Variable Gain Photoreceiver – Fast Optical Power Meter

Related OE-200 Models

@ 850 nm

- OE-200-SI-FST      Si-PIN, Ø 1.2 mm, 320 - 1060 nm  
free space input, 1.035"-40 threaded flange
- OE-200-SI-FS      Si-PIN, Ø 1.2 mm, 320 - 1060 nm  
free space input, 25 mm dia. unthreaded flange
- OE-200-SI-FC      Si-PIN, Ø 1.2 mm, 320 - 1060 nm  
FC fiber connector (fix/permanent)

@ 1310 nm

- OE-200-IN1-FST    InGaAs-PIN, Ø 300 µm, 900 - 1700 nm  
free space input, 1.035"-40 threaded flange
- OE-200-IN1-FS    InGaAs-PIN, Ø 300 µm, 900 - 1700 nm  
free space input, 25 mm dia. unthreaded flange
- OE-200-IN1-FC    InGaAs-PIN, integrated ball lens, 900 - 1700 nm  
FC fiber connector (fix/permanent)

@ 1550 nm

- OE-200-IN2-FST    InGaAs-PIN, Ø 300 µm, 900 - 1700 nm  
free space input, 1.035"-40 threaded flange
- OE-200-IN2-FS    InGaAs-PIN, Ø 300 µm, 900 - 1700 nm  
free space input, 25 mm dia. unthreaded flange
- OE-200-IN2-FC    InGaAs-PIN, integrated ball lens, 900 - 1700 nm  
FC fiber connector (fix/permanent)

OE-200-S      customized versions available on request

Available Accessories

PRA-FSMA  
PRA-FC



fiber-adapter with external  
1.035"-40 thread

PRA-PAP



post adapter plate,  
easy to mount on  
FEMTO photoreceiver series OE,  
FWPR, PWPR, HCA-S and LCA-S

PS-15




power supply,  
input: 100 - 240 VAC,  
output: ±15 VDC

LUCI-10



compact digital I/O interface for USB remote control,  
supports opto-isolation of amplifier signal path from PC  
USB port, 16 digital outputs, 3 opto-isolated digital inputs,  
bus-powered operation

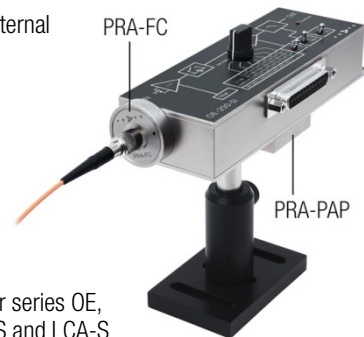

## Variable Gain Photoreceiver – Fast Optical Power Meter

Specifications	<p>Test conditions <math>V_s = \pm 15\text{ V}</math>, <math>T_A = 25\text{ }^\circ\text{C}</math>, output load impedance <math>1\text{ M}\Omega</math></p>
Gain	<p>Conversion gain <math>1 \times 10^3 \dots 1 \times 10^{11}\text{ V/W}</math> (@ 850 nm, output load <math>\geq 100\text{ k}\Omega</math>)</p> <p>Gain accuracy <math>\pm 1\%</math> electrical, between settings</p> <p>Conversion gain accuracy</p> <p style="margin-left: 20px;">OE-200-UV-FST/FS (@ <math>P_{\text{OPT}} \leq 2\text{ mW}</math>, 850 nm) free space <math>\pm 15\%</math> nominal</p> <p style="margin-left: 20px;">OE-200-SI-FST (@ <math>P_{\text{OPT}} \leq 2\text{ mW}</math>, 1550 nm) with fiber adapter (PRA series) <math>\pm 15\%</math> nominal</p> <p style="margin-left: 20px;">OE-200-UV-FC (@ <math>P_{\text{OPT}} \leq 1\text{ mW}</math>, 850 nm) fixed fiber input connector <math>\pm 5\%</math> guaranteed by factory calibration*</p> <p>* Factory verified with MM 50/125, FC/APC, NA 0.22 (when using FC/PC fiber connector, coupling efficiency may differ slightly.) Coupling efficiency depends on fiber type.</p> <p>Gain drift see table below</p>
Frequency Response	<p>Lower cut-off frequency DC / 1 Hz, switchable</p> <p>Upper cut-off frequency (-3dB) up to 500 kHz (see table below), switchable to 10 Hz</p>
Detector	<p>Detector type Si-PIN photodiode</p> <p>Active area <math>1.1 \times 1.1\text{ mm}^2</math></p> <p>Spectral range 190 - 1000 nm</p> <p>Sensitivity</p> <p style="margin-left: 20px;"><math>0.29\text{ A/W}</math> (@ 850 nm)</p> <p style="margin-left: 20px;"><math>0.36\text{ A/W}</math> (@ 700 nm)</p>
Input	<p>Input offset current (dark current) 2 pA typ.</p> <p>Input offset drift see table below</p> <p>Input offset compensation range <math>\pm 600\text{ pA}</math>, adjustable by offset potentiometer or <math>\pm 400\text{ pA}</math>, adjustable by external control voltage</p> <p>Optical CW saturation power see table below</p> <p>Noise equivalent power (NEP) see table below</p>

## Variable Gain Photoreceiver – Fast Optical Power Meter

Specifications (continued)

Performance Depending  
on Gain Setting

Gain setting (low noise) (V/W)**	10 <sup>3</sup>	10 <sup>4</sup>	10 <sup>5</sup>	10 <sup>6</sup>	10 <sup>7</sup>	10 <sup>8</sup>	10 <sup>9</sup>
Upper cut-off frequency (-3 dB)	500 kHz	500 kHz	400 kHz	200 kHz	50 kHz	7 kHz	1.1 kHz
Rise/fall time (10 % - 90 %)	700 ns	700 ns	900 ns	1.8 μs	7 μs	50 μs	300 μs
NEP (√Hz)**	60 pW	7.3 pW	1.5 pW	450 fW	150 fW	48 fW	17 fW
Measured at	10 kHz	10 kHz	10 kHz	1 kHz	1 kHz	100 Hz	100 Hz
Integr. input noise (RMS)***	63 nW	9 nW	2.8 nW	1 nW	320 pW	46 pW	6.2 pW
Input offset drift (°C)**	100 nW	10 nW	1 nW	85 pW	8.5 pW	1.3 pW	1 pW
Gain drift (°C)	0.008%	0.008%	0.008%	0.01%	0.01%	0.01%	0.02%
Optical CW saturation power**	2 mW	1 mW	0.1 mW	10 μW	1 μW	0.1 μW	10 nW

Gain setting (high speed) (V/W)**	10 <sup>5</sup>	10 <sup>6</sup>	10 <sup>7</sup>	10 <sup>8</sup>	10 <sup>9</sup>	10 <sup>10</sup>	10 <sup>11</sup>
Upper cut-off frequency (-3 dB)	500 kHz	500 kHz	400 kHz	200 kHz	50 kHz	7 kHz	1.1 kHz
Rise/fall time (10 % - 90 %)	700 ns	700 ns	900 ns	1.8 μs	7 μs	50 μs	300 μs
NEP (√Hz)**	48 pW	6.6 pW	1.5 pW	450 fW	150 fW	48 fW	17 fW
Measured at	10 kHz	10 kHz	10 kHz	1 kHz	1 kHz	100 Hz	100 Hz
Integr. input noise (RMS)***	41 nW	6.8 nW	2.5 nW	920 pW	300 pW	43 pW	6.1 pW
Input offset drift (°C)**	100 nW	10 nW	1 nW	85 pW	8.5 pW	1.3 pW	1 pW
Gain drift (°C)	0.008%	0.008%	0.008%	0.01%	0.01%	0.01%	0.02%
Optical CW saturation power**	0.1 mW	10 μW	1 μW	0.1 μW	10 nW	1 nW	0.1 nW

\*\* referred to 850 nm

\*\*\* The integrated input noise is measured with a shaded input in the full bandwidth ("FBW") setting (referred to 850 nm).

The input referred peak-peak noise can be calculated from the RMS noise as follows:

$$P_{\text{Input noise peak-to-peak}} = P_{\text{Input noise RMS}} \times 6$$

The output noise is given by:

$$U_{\text{Output noise RMS}} = P_{\text{Input noise RMS}} \times \text{gain}$$

$$U_{\text{Output noise peak-to-peak}} = U_{\text{Output noise RMS}} \times 6 = P_{\text{Input noise RMS}} \times \text{gain} \times 6$$

The integrated noise will be reduced considerably by setting the low pass filter to "10 Hz" instead of "FBW". This is especially useful for continuous wave (CW) measurements.

Output

Output voltage range	±10 V (@ ≥100 kΩ output load)
Max. output current	±30 mA (short-circuit proof)
Output impedance	50 Ω (terminate with ≥100 kΩ)

Indicator LED

Function	overload
----------	----------

Digital Control

Control input voltage range	LOW bit: -0.8 ... +1.2 V, HIGH bit: +2.3 ... +12 V
Control input current	0 mA @ 0 V, 1.5 mA @ +5 V, 4.5 mA @ +12 V
Overload output	nonactive: <0.4 V, @ 0 ... -1 mA active: typ. 5 ... 5.1 V @ 0 ... 2 mA

Ext. Offset Control

Control voltage range	±10 V
Offset control input impedance	20 kΩ
Conversion factor	40 pA/V

Power Supply

Supply voltage	±15 V (±14.75 ... ±16.5 V)
Supply current	+110/-80 mA (depends on operating conditions, recommended power supply capability min. ±200 mA)
Stabilized power supply output	±12 V, max. 50 mA, +5 V, max. 30 mA

Case

Weight	360 g (0.79 lb)
Material	AlMg4.5Mn, nickel-plated

Temperature Range

Storage temperature	-40 ... +80 °C
Operating temperature	0 ... +60 °C

## Variable Gain Photoreceiver – Fast Optical Power Meter

<p>Absolute Maximum Ratings</p>	<p>Optical input power (CW)            20 mW                  Digital control input voltage        –5 V/+16 V relative to digital ground DGND (pin 9)                  Analog control input voltage       ±15 V relative to analog ground AGND (pin 3)                  Power supply voltage                ±20 V</p>
<p>Connectors</p>	<p>Input</p> <p style="margin-left: 20px;">OE-200-UV-FST      1.035" -40 threaded flange for free space applications</p> <p style="margin-left: 20px;">OE-200-UV-FS        25 mm unthreaded flange for free space applications</p> <p style="margin-left: 20px;">OE-200-UV-FC        FC fiber optic connector</p> <p>Output</p> <p style="margin-left: 20px;">BNC jack (female)</p> <p>Power supply</p> <p style="margin-left: 20px;">Lemo® series 1S, 3-pin fixed socket (mating plug type: FFA.1S.303.CLAC52)                  Pin 1:    +15 V                  Pin 2:    –15 V                  Pin 3:    GND</p> <div style="text-align: center; margin: 10px 0;"> </div> <p>Control port</p> <p style="margin-left: 20px;">Sub-D 25-pin, female, qual. class 2                  Pin 1:    +12 V (stabilized power supply output)                  Pin 2:    –12 V (stabilized power supply output)                  Pin 3:    AGND (analog ground for pins 1 - 8)                  Pin 4:    +5 V (stabilized power supply output)                  Pin 5:    overload output: HIGH = overload (referred to pin 3)                  Pin 6:    signal output (connected to BNC)                  Pin 7:    NC                  Pin 8:    input offset control voltage                  Pin 9:    DGND (ground for digital control pins 10 - 14)                  Pin 10:   digital control input: gain, LSB                  Pin 11:   digital control input: gain                  Pin 12:   digital control input: gain, MSB                  Pin 13:   digital control input: AC/DC                  Pin 14:   digital control input: high speed / low noise                  Pin 15 - 25: NC</p>
<p>Scope of Delivery</p>	<p>OE-200-UV, internally threaded coupler ring (FST version only), Lemo® 3-pin connector, datasheet, transport package</p>

## Variable Gain Photoreceiver – Fast Optical Power Meter

Remote Control Operation

General

Remote control input bits are opto-isolated and connected by a logical OR function to the local switch settings. For remote control set the corresponding local switches to “Remote”, “AC” and “H” and select the desired setting via a bit code at the corresponding digital inputs. Mixed operation, e.g. local AC/DC setting and remote controlled gain setting, is also possible. The switch setting “FBW / 10 Hz” of the low pass signal filter is not remote controllable.

Gain setting

Low noise Gain (V/W) Pin 14=HIGH	High speed Gain (V/W) Pin 14=LOW	Pin 12 MSB	Pin 11	Pin 10 LSB
$10^3$	$10^5$	LOW	LOW	LOW
$10^4$	$10^6$	LOW	LOW	HIGH
$10^5$	$10^7$	LOW	HIGH	LOW
$10^6$	$10^8$	LOW	HIGH	HIGH
$10^7$	$10^9$	HIGH	LOW	LOW
$10^8$	$10^{10}$	HIGH	LOW	HIGH
$10^9$	$10^{11}$	HIGH	HIGH	LOW

Gain settling time


<150 ms

AC/DC setting

Coupling	Pin 13
AC	LOW
DC	HIGH

Conversion Gain


Normalized Conversion Gain

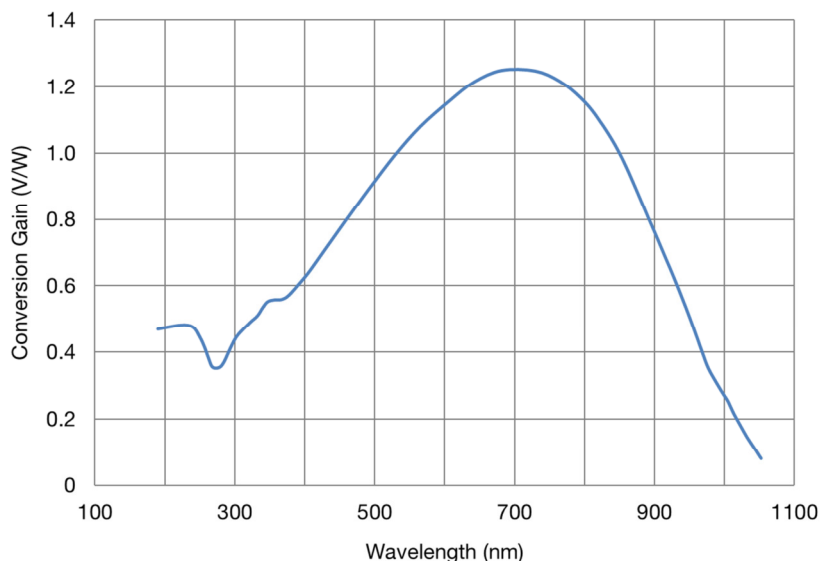


# Variable Gain Photoreceiver – Fast Optical Power Meter

Dimensions

OE-200-UV-FST (1.035"-40 threaded free space input):

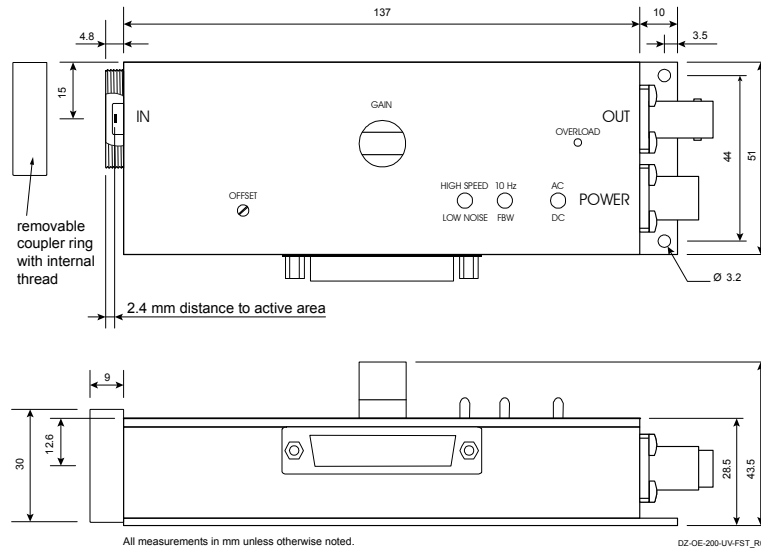

OE-200-UV-FS (25 mm dia. unthreaded free space input):

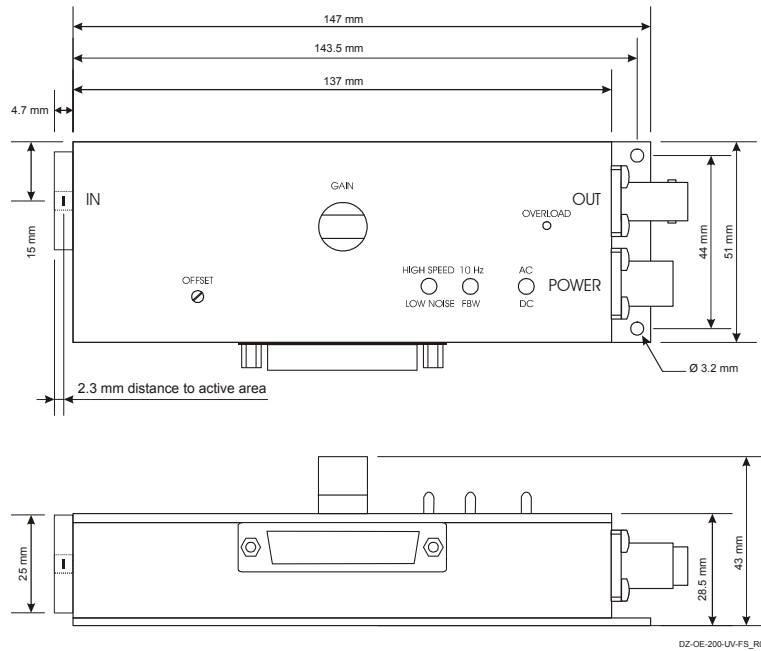




# Variable Gain Photoreceiver – Fast Optical Power Meter

Dimensions (continued)

OE-200-UV-FC (FC fiber optic input):

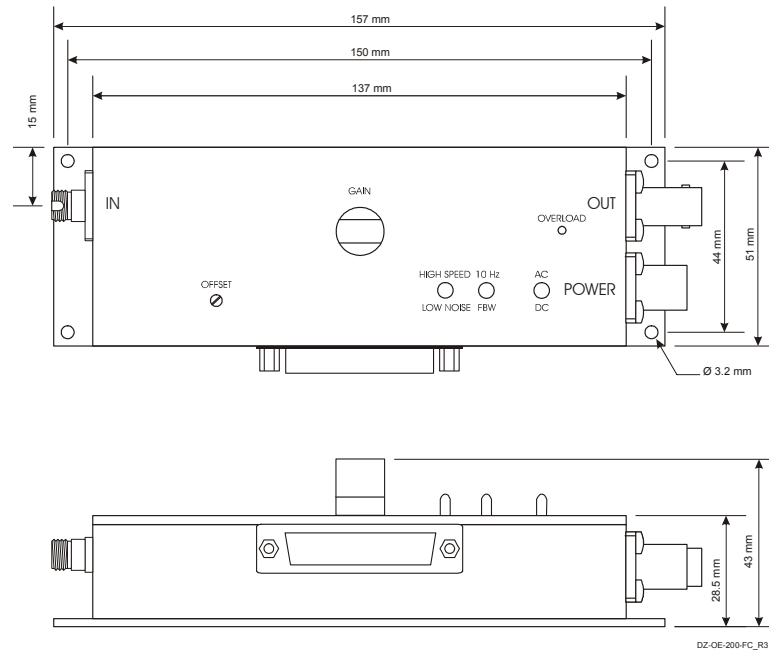

Specifications are subject to change without notice. Information provided herein is believed to be accurate and reliable. However, no responsibility is assumed by FEMTO Messtechnik GmbH for its use, nor for any infringement of patents or other rights of third parties which may result from its use. No license is granted by implication or otherwise under any patent or patent rights of FEMTO Messtechnik GmbH. Product names mentioned may also be trademarks used here for identification purposes only.

© by FEMTO Messtechnik GmbH · Printed in Germany